

Sequencing using the ABI 377

1. Pour gel.
2. Damp paper towel (~0.5 hr)
3. Wash plates and allow to dry.
4. Put plates in mounting rack.
5. Add bottom buffer tray.
6. Put on plates.
7. Do plate check.
8. Fill bottom buffer tray.
9. Add heating plate.
10. Do pre-run.
11. Prepare samples
12. Load Comb.
13. Fill well with ficoll.
14. Put in comb.
15. Put on upper buffer chamber and fill.
16. Run for 1 minute.
17. Remove comb.
18. Rinse out ficoll.
19. Put on upper buffer chamber lid.
20. Resume run.
21. Check status and log.
22. Clean comb.
23. Clean work area.